



IX81 specifications

Microscope body	Revolving nosepiece		Sextuple motorized with objective lens retraction in PC mode, simple waterproof mechanism incorporated
	Focus		9mm stroke (from stage surface, 7mm upward and 2mm downward), fine/coarse switchable focusing knobs (minimum graduation: 0.1µm), objective lens escape/return buttons and return to memory position buttons (each side of microscope frame)
	Primary image port		Lower port (standard left side port: S1F 100% or S8F 80%, or optional lower Back port selectable, 2-step light path selection), Upper port when built-in magnification changer 1x/1.6x is replaced (optional right side port or upper back port selectable, 2-step light path selection), Bottom port (option)
	Frontal operation		Light path selector button, Transmitted light intensity control buttons and light ON/OFF switch button, Fine/Coarse focus selector button, TTL Pulse control switch (auxiliary) buttons
Transmitted light illuminator	100W transmitted light illumination pillar	IX2-ILL100	Pillar tilt mechanism (30° inclination angle, with shock decreasing mechanism), Condenser holder (with 50mm stroke, swing-in/out mechanism), Field iris diaphragm adjustable, 4 filter holders (ø45mm, t=6mm or less)
	External power supply unit	IX2-UCB	Auto voltage selector (100V / 200V), RS232C interface for PC operation, IX2-BSW driver software
Observation tube	Tilting binocular tube	U-TBI90	35-85° continuous angle adjustable (eyepoint height range: 406mm-471mm), interpupillary distance adjustable between 50-76mm, diopter adjustment function, erect image, F.N. 22
	Binocular tube	U-BI90CT	Built-in focusing telescope, interpupillary distance adjustable 50-76mm, diopter adjustment function, F.N. 22
		U-BI90	Interpupillary distance adjustable 50-76mm, diopter adjustment function, F.N. 22
Trinocular tube	TR30H+IX-ATU	3 step optical path selectable (observation: straight port = 100:0, 20:80, 0:100), interpupillary distance adjustable 50-76mm, diopter adjustment function, F.N. 22	
Stage	Cross stage with flexible right handle	IX2-SFR	50mm(X) x 50mm(Y) stroke, stage insert plate exchangeable (ø110mm)
	Cross stage with short left handle	IX-SVL-2	50mm(X) x 43mm(Y) stroke, stage insert plate exchangeable (ø110mm)
	Plain stage	IX2-SP	232mm(X) x 240mm(Y) stage size, stage insert plate exchangeable (ø110mm)
		IX-MVR	Mechanical stage to be used with IX2-SP, 130mm(X) x 85mm(Y) stroke
	Narrow plain stage	IX2-KSP	160mm(X) x 240mm(Y) stage size, stage insert plate exchangeable (ø110mm)
		CK40-MVR	Mechanical stage to be used with IX2-KSP, 120mm(X) x 78mm(Y) stroke
Gliding stage	IX2-GS	Upper circular stage 360° rotatable, 20mm(X/Y) travel	
Condenser	Motorized long working distance universal	IX2-LWUCDA	Motorized turret with 6 position slots for optical devices (3 positions each for ø30mm and ø38mm), aperture iris diaphragm adjustable, N.A. 0.55 / W.D. 27mm
	Long working distance Relief Contrast	IX2-MLWCD	4 positions for optical devices (for ø50mm, Relief Contrast optical devices rotatable), aperture iris diaphragm adjustable, N.A. 0.5 / W.D. 45mm
	Ultra long working distance	IX-ULWCD	4 positions for optical devices (for ø29mm), aperture iris diaphragm adjustable, N.A. 0.3 / W.D. 73mm
	Water immersion DIC	IX2-DICD+IX2-TLW	Single position for optical device (include two optical device holders), 40 injection pipette or electrode insertion angle, aperture iris diaphragm adjustable, N.A. 0.9 / W.D. 3.7mm
Eyepiece		WH10x	High eyepoint, F.N. 22
		WH10x-H	High eyepoint, diopter adjustment function, F.N. 22
Reflected light fluorescence unit	Fluorescence illuminator	IX2-RFAL	L-shaped design with exchangeable F.S. and A.S. modules, two filter holder sliders (2 positions, ø32mm, t=6mm or less)
		IX2-RFA	Straight design with field iris diaphragm, filter holder slider (2 positions, ø32mm, t=6mm or less)
	Fluorescence cube turret	IX2-RFACA	Motorized turret with 6 positions, built-in shutter
	Light source		100W HBO lamp housing and HBO transformer, or 75WXBO lamp housing and XBO transformer


IX81 dimensions

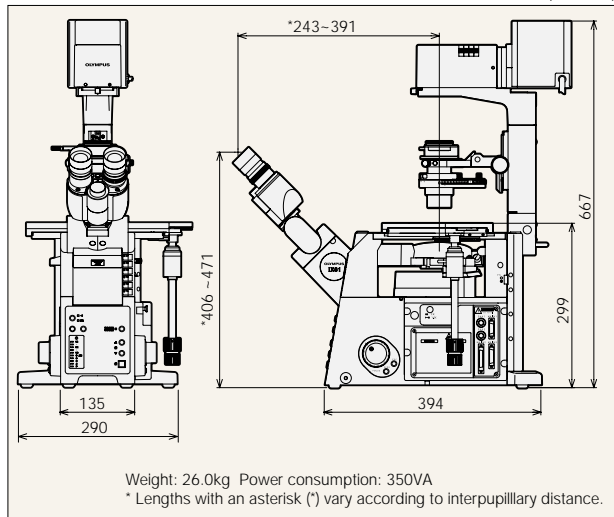


IX2-UCB dimensions

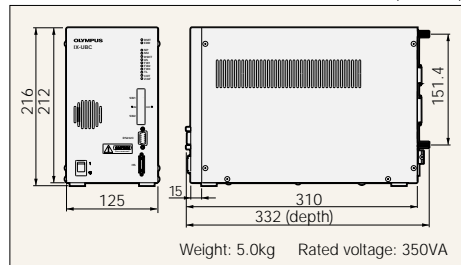

U-FH dimensions

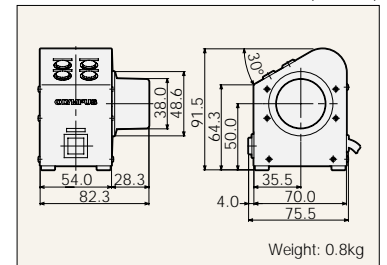

U-HSTR2 dimensions

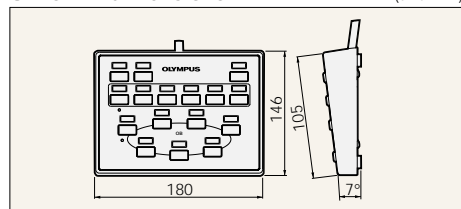

Photos courtesy of : Dr. Hiraoka, Cell Biology Group Leader, Kansai Advanced Research Center, Communications Research Laboratory (P1-2 time-lapse)
Pr. Yasuhiro Kamei, Pr. Shunsuke Yuba, Institute for Molecular and Cellular Biology, Osaka University
(P8 below, coexpression of EGFP and DsRed in a Zebrafish embryo, extended focus image of 10µmX28 slice)


Specifications are subject to change without any obligation on the part of the manufacturer.



OLYMPUS

OLYMPUS OPTICAL CO., LTD.
San-Ei building, 22-2, Nishi Shinjuku 1-chome, Shinjuku-ku, Tokyo, Japan
OLYMPUS OPTICAL CO. (EUROPA) GMBH.
Postfach 10 49 08, 20034, Hamburg, Germany
OLYMPUS AMERICA INC.
2 Corporate Center Drive, Melville, NY 11747-3157, U.S.A.
OLYMPUS SINGAPORE PTE LTD.
491B River Valley Road, #12-01/04 Valley Point Office Tower, Singapore 248373
OLYMPUS OPTICAL CO. (U.K.) LTD.
2-8 Honduras Street, London EC1Y 0TX, United Kingdom.
OLYMPUS AUSTRALIA PTY. LTD.
104 Ferntree Gully Road, Oakleigh, Victoria, 3166, Australia

OLYMPUS BUSINESS AREAS

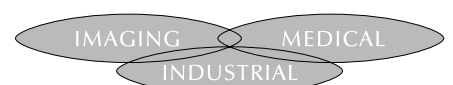

www.olympus.com

Printed in Japan M3502E-0302B

OLYMPUS[®]


MOTORIZED
INVERTED SYSTEM MICROSCOPE


IX81
IX2 SERIES

Automated Platform for
Live Cell Imaging

UIS
UNIVERSAL
INFINITY SYSTEM



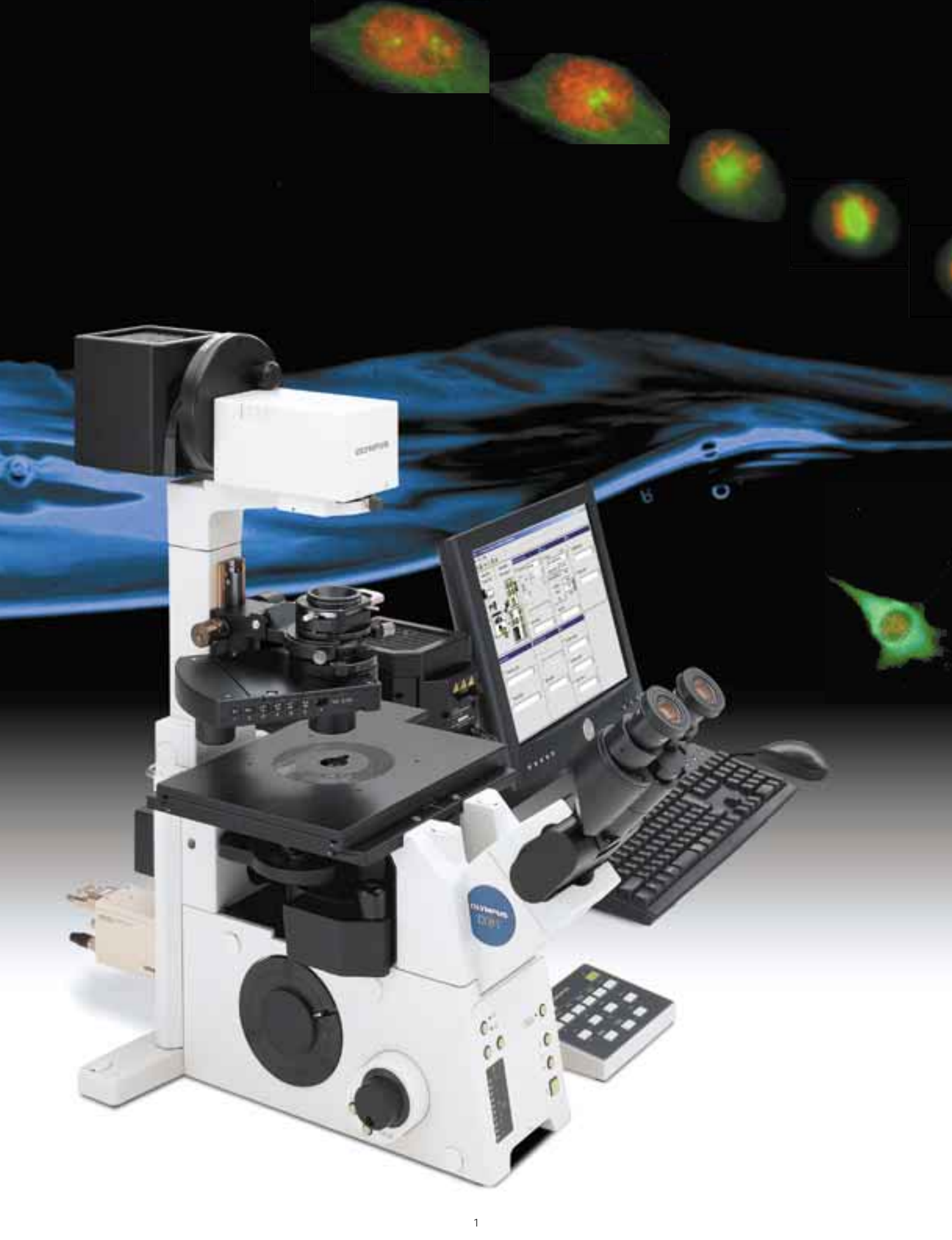

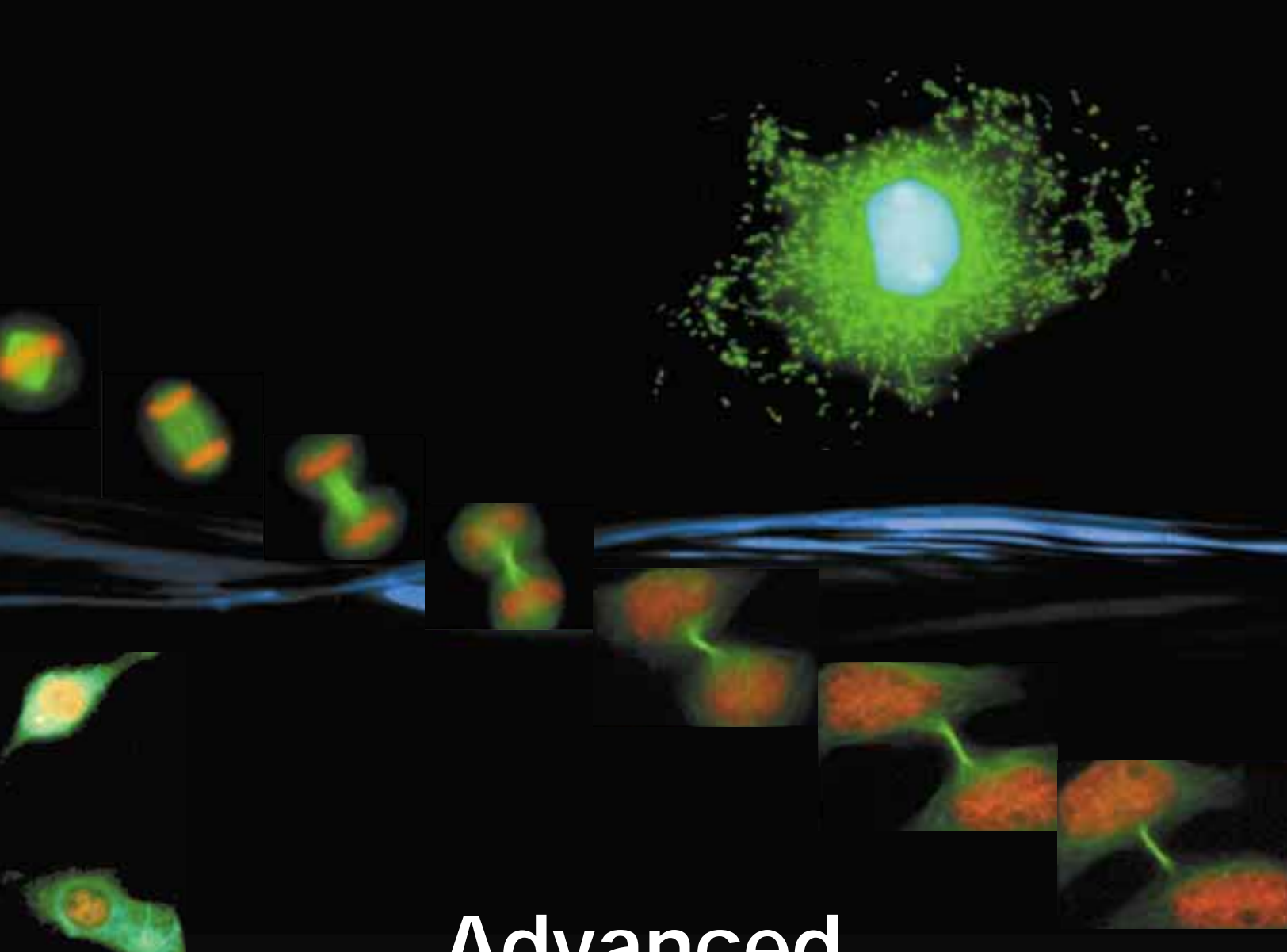

Advanced, fully motorized system for superior performance.

The IX81 is Olympus' most advanced motorized microscope, providing high standards of observation, measurement and functional operation. A V-shaped optical path, improved fluorescence illuminators and further expanded UIS optics provide excellent performance for research applications. A modular system of motorized accessories allows for complete customization to individual research needs. Multi-dimensional analyses can be performed automatically or even remotely via computer control. Additionally, the IX81 features numerous input and output ports allowing the use of several light sources and detectors. The IX81 is the optimum imaging platform for research and automation.

**Note:*

UIS optics: The original Olympus infinity-corrected optical system. This system expresses the superb Olympus optical technology, provides high resolution, high contrast images and the flexibility to meet current and future applications.

A flexible motorized platform for automated system solutions.

A selection of independent motorized units for convenient operation.

All major microscope functions can be motorized, including focus, illumination, objectives and optical path selection. The motorized units for the IX81 can be operated via a remote handset or buttons on the microscope body. All controls are close to the operator's hand, maximizing user comfort while minimizing needed bench space.

Computer control for routine and research applications.

All motorized functions are also controllable via computer. Imaging software can be used to facilitate microscope operations ranging from routine to complex. Buttons on the microscope body and remote handset can be programmed for custom operation.



Motorized universal condenser/IX2-LWUCDA

6 turret positions allow brightfield, phase contrast and DIC observations. Also accepts 32mm filters.



Handset/U-HSTR2

A remote handset controls all motorized functions via a convenient and programmable interface.



Focus handle/U-FH

The remote focus handle duplicates the feel and function of the microscope's focus knobs. Additional controls include fine/coarse focus selection, lamp on/off, shutter open/close, and camera vs. visual observation.




System controller/IX2-UCB

All motorized units are powered by this external system controller. Included is an RS232C connection for PC commands and expansion slots for future system upgrades.



Microscope front panel

Easy to use buttons allow selection of light path, focus fine/coarse switchover, light intensity and lamp on/off control. Two auxiliary buttons can be custom programmed. Includes LED lamp intensity meter.

- 
- Light path selector
 - Auxiliary buttons
 - Fine/Coarse
 - Transmitted light intensity control
 - Transmitted light ON/OFF



Motorized shutter/IX2-SHA

Can be mounted in both transmitted and reflected light paths.



**Motorized filter wheel/
U-FWR and U-FWO**

6 filter positions
(32mm or 25mm diameter).



Motorized sextuple revolving nosepiece

Included with microscope frame.

**Motorized fluorescence cube turret/
IX2-RFACA**

6 fluorescence filter cube positions.



Internal motorized focus drive

0.01 μm minimum step size.

**Objective escape and
zero-return buttons**

Moves objective to lower focus limit.
Allows setting of default focus position.

* Equipped on each side of microscope frame.



**Motorized bottom port unit with C-mount/
IX2-TVRAC**




Escape button

Return to memory position button

Unique two-tiered multi-port design offers input/output system flexibility and application support

The IX frame begins with 6 input/output ports for a wide variety of light sources or cameras. The unique Olympus design allows up to four ports to have simultaneous access to the primary image, without relay lenses, providing cameras with the highest quality image, free of aberrations. Two-tiered multi-port design maximizes flexibility while maintaining a 299mm low stage height. The compact and rigid frame requires a minimum amount of bench space, easily accommodates external equipment and remains stable during extended observations. Over 100mm of free space is available at the left side port between the frame and the primary image, easily allowing for adaptation of instrumentation for individual research requirements.

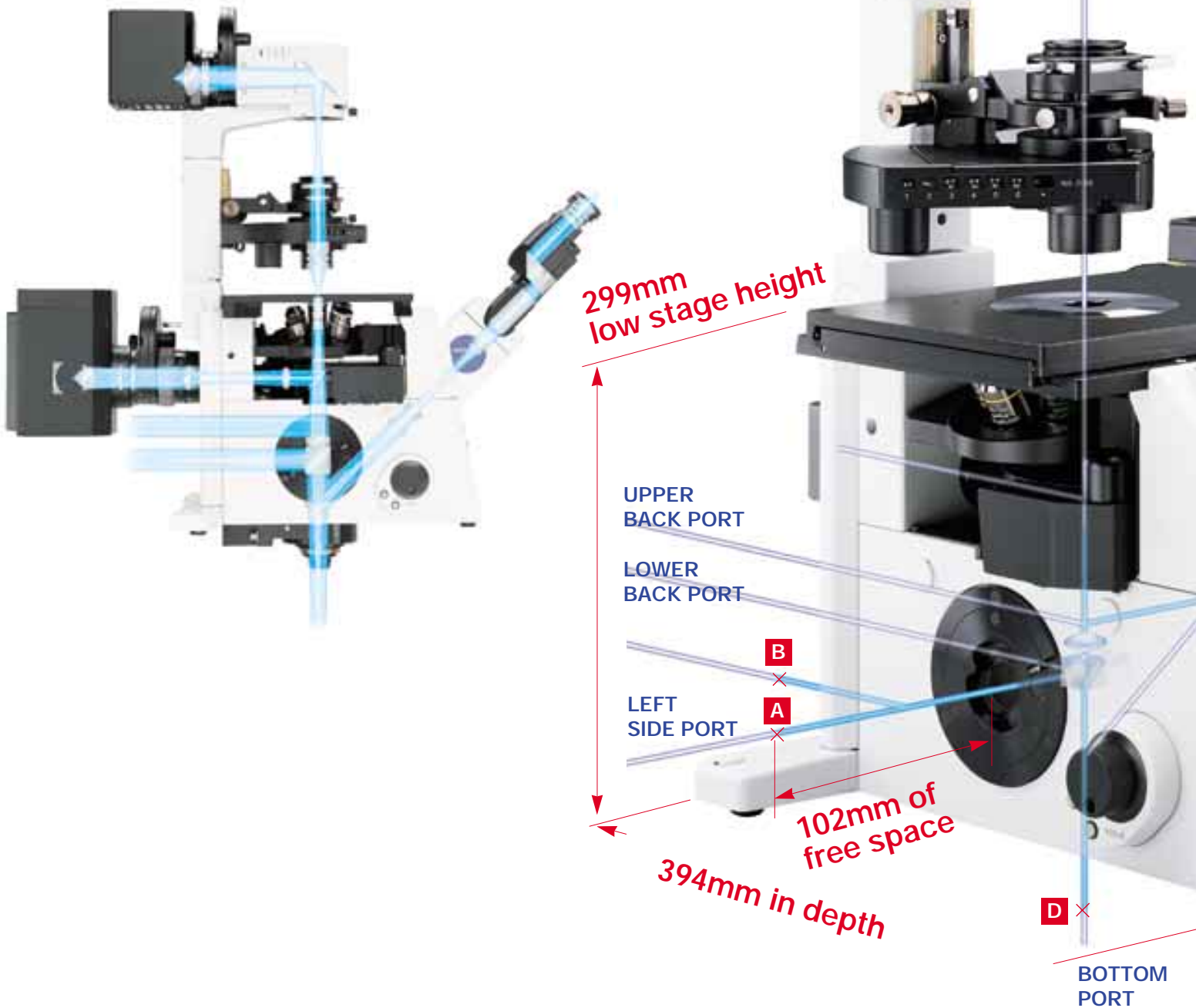

A B C D

Primary images are obtainable from all 4 ports simultaneously



LEFT SIDE PORT (Primary Image)

A B

For film or CCD camera systems. Double port tube allows the attachment of two cameras (both primary images). Also accepts Olympus confocal scan units.



U-DPCAD double port tube

RIGHT SIDE PORT (Primary Image)

C

Used for connecting an additional light source or CCD camera.



IX2-RSPC
Right side port attachment with C-mount

BOTTOM PORT (Primary Image)

D

Well suited for extremely sensitive low light level imaging.



IX2-TVR bottom port adapter
+ CCD camera

UPPER BACK PORT (Primary Image)

Used for connecting an additional light source or CCD camera.

LOWER BACK PORT (Primary Image)

For the attachment of large bodied items such as cooled CCD cameras or confocal scan units.



BINOCULAR PORT

CCD camera mount via side port intermediate tube.



GX-SPU
side port
intermediate
tube

High quality images meet the demands of advanced analytical applications.

New fluorescence accessories for improved sensitivity

Improvements in the performance of the UIS optical system have resulted in sharper, clearer images. Two reflected light fluorescence illuminators are now available.

- An L-shaped illuminator includes centerable aperture and field stops with convenient controls close to the user. L-shape also permits the mounting of a large bodied camera to the back port.
- A straight illuminator designed for maximum intensity provides the brightest possible images.



A variety of fluorescence filter cubes are available. Up to 6 cubes can be installed in the motorized turret, with easy selection via the remote handset. An optional DIC filter cube incorporates an analyzer and allows automatic switching between DIC observation and fluorescence.

Rectangular field stop/IX2-RFSS

Pinhole field stop/IX2-RFSPOT

Flexible field stop options IX2-RFSS rectangular field stop module or IX2-RFSPOT pinhole field stop module can be mounted into the



L-shaped illuminator for improved illumination control during fluorescence imaging or photobleaching experiments.

(Available autumn 2002)


GFP, CFP, YFP Filter Cubes

Fluorescence performance has been improved with high efficiency coatings minimizing signal crossover. Filter sets are optimized for



specific fluorochromes such as GFP, CFP and YFP. All Olympus filter cube housings incorporate anti-reflective materials that minimize stray light .

(Available autumn 2002)

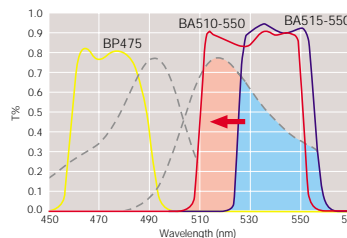

The new IBA cube has a capability to maximize deflection of GFP emission.

Long working distance objectives/LUCPLFL40x

Accommodating vessel thicknesses from 0-2mm with a correction collar mechanism that maintains objective focus during adjustment.



Correcting for different vessels is fast and simple. Lens design for 340nm transmission allows these objectives to be used for UV fluorescence applications such as FURA-2.

High quality images of living cells

The IX2-LWUCDA motorized universal condenser enables standard brightfield, phase contrast and DIC observations. Three series of DIC prisms are available for optimum contrast enhancement with a wide variety of specimens. Standard prisms are best for most



observations, while High Contrast (HC) prisms provide maximum contrast for thin specimens, and High Resolution (HR) prisms provide maximum detail with thicker specimens.

Optimum images for Microinjection


Relief Contrast Condenser provides long working distance (W.D. 45mm) for injector access and a high numerical aperture (N.A. 0.5) for excellent images with plastic dishes. With a long working distance and high N.A. Phase contrast and DIC are also possible.



Unique Relief Contrast objectives feature rotatable modulators for consistent shading for all magnifications.

Computer controlled multi-color and time-lapse applications


Motorized precise internal Z-drive, objective nosepiece and accessories allow multi-dimensional, multi-spectral imaging via a computer. A thermally and structurally stable frame design combine to create an ideal system for long term time-lapse and deconvolution observations.

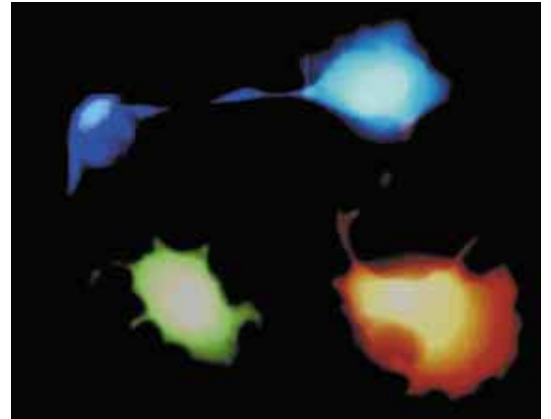


Fully automatic confocal system to meet a wide range of needs


The Fluoview confocal scan unit can be attached to the back port, leaving ample space for manipulators and electrophysiological equipment. Combining the IX81 with the Fluoview FV500 provides a fully capable, multi-channel, motorized system. Both confocal and fluorescence observation equipment can be attached simultaneously, and port changing can be performed automatically.

*Special microscope body available for Fluoview.

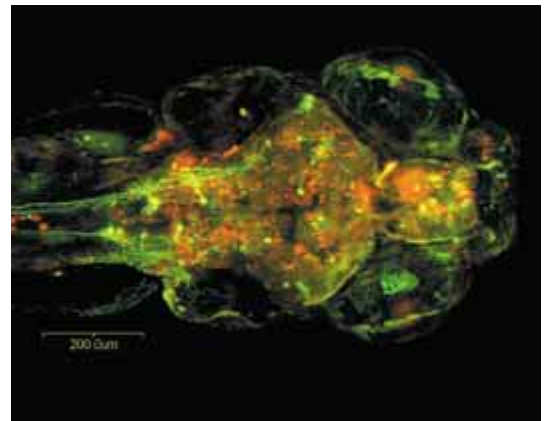

IX81 SYSTEM DIAGRAM

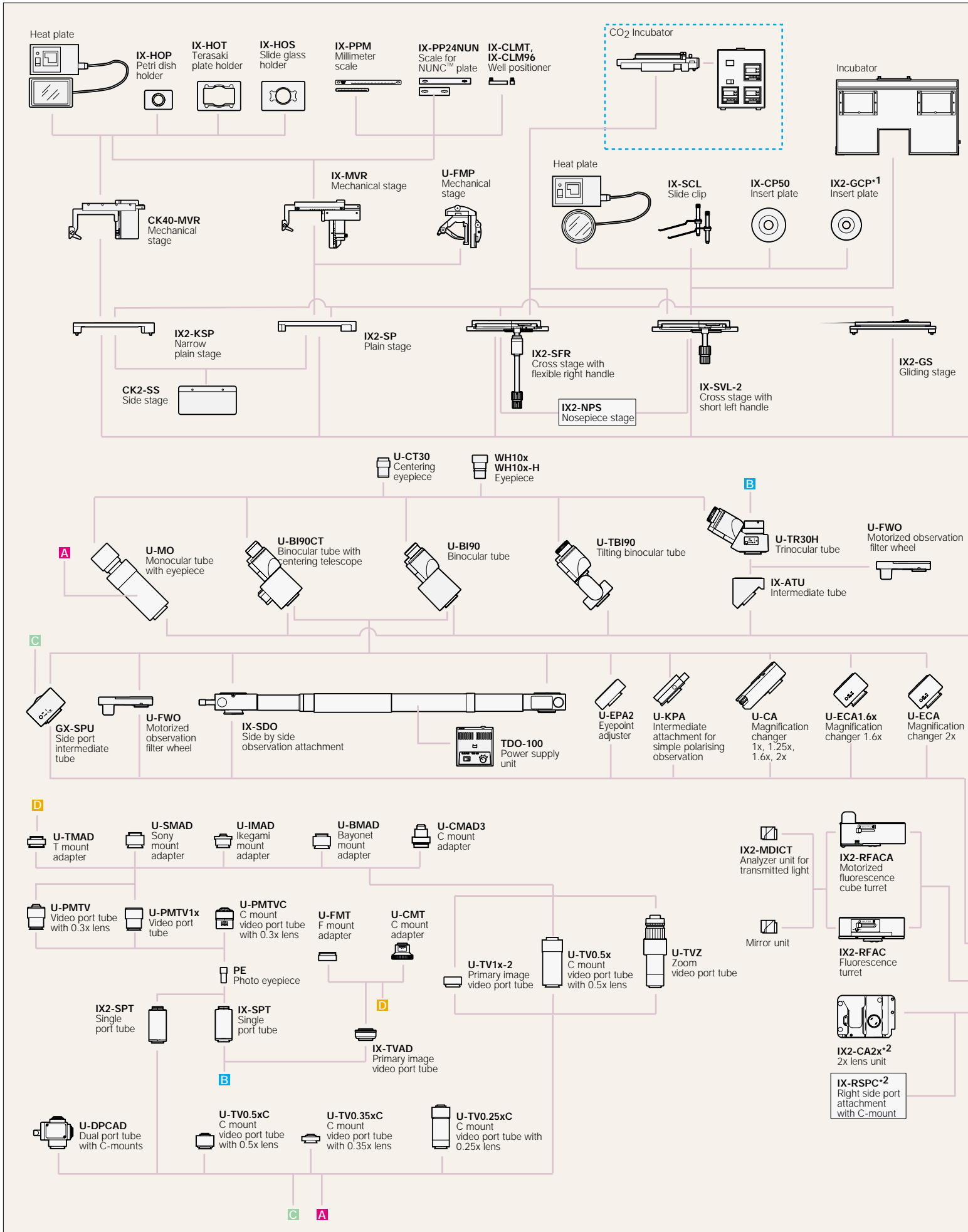

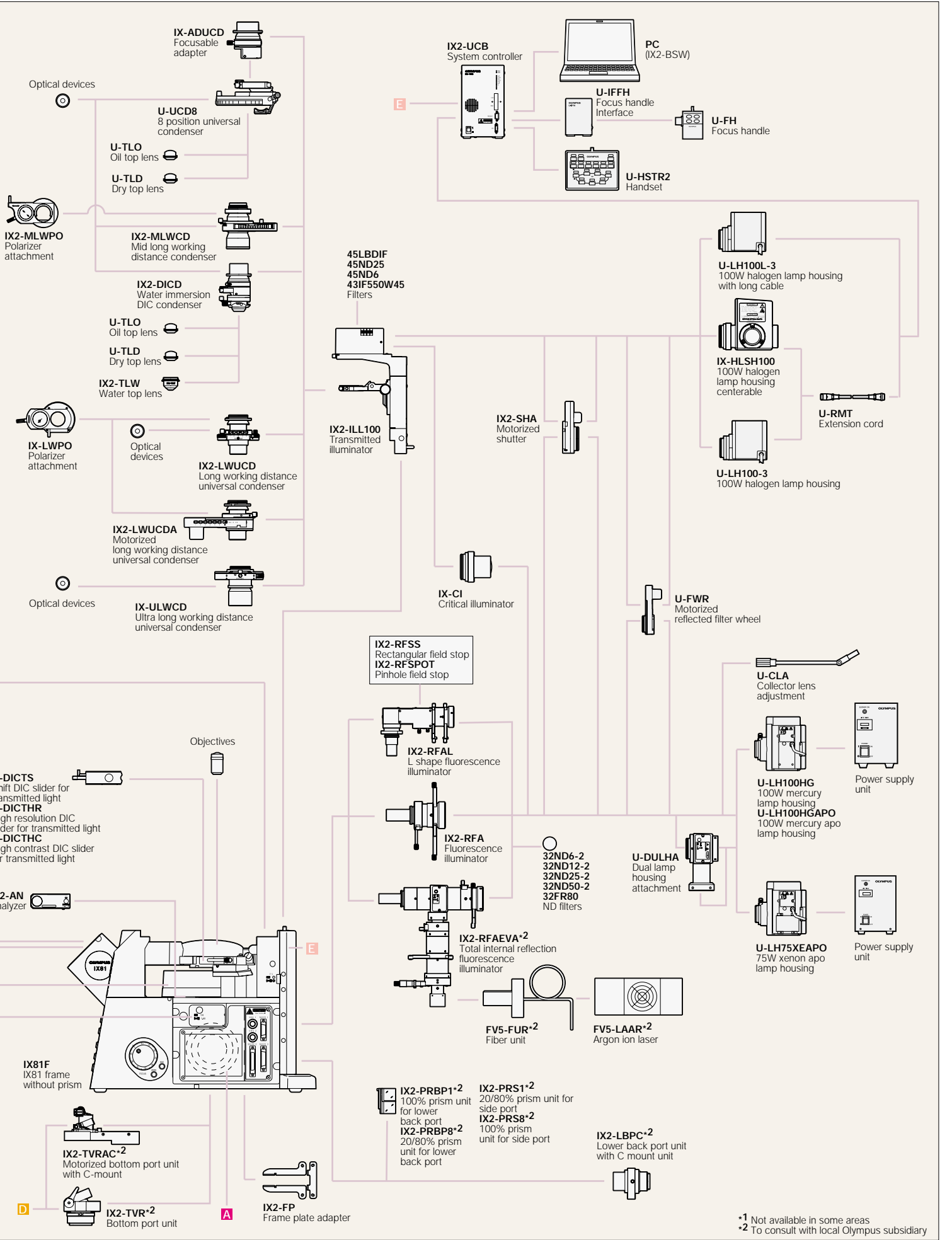

IX-ADUCD
Focusable adapter

U-UCD8
8 position universal condenser

U-TLO
Oil top lens

U-TLD
Dry top lens

IX2-MLWPO
Polarizer attachment

IX2-MLWCD
Mid long working distance condenser

IX2-DICD
Water immersion DIC condenser

U-TLO
Oil top lens

U-TLD
Dry top lens

IX2-TLW
Water top lens

IX-LWPO
Polarizer attachment

IX2-LWUCD
Long working distance universal condenser

IX2-LWUCDA
Motorized long working distance universal condenser

IX2-ULWCD
Ultra long working distance universal condenser

U-DICTS
Shift DIC slider for transmitted light

U-DICTHR
High resolution DIC slider for transmitted light

U-DICTHC
High contrast DIC slider for transmitted light

IX2-AN
Analyzer

IX81F
IX81 frame without prism

IX2-TVRAC*2
Motorized bottom port unit with C-mount

IX2-TVR*2
Bottom port unit

IX2-FP
Frame plate adapter

IX2-UCB
System controller

PC
(IX2-BSW)

U-IFFH
Focus handle Interface

U-FH
Focus handle

U-HSTR2
Handset

45LBDIF
45ND25
45ND6
43IF550W45
Filters

IX2-ILL100
Transmitted illuminator

IX2-SHA
Motorized shutter

U-LH100L-3
100W halogen lamp housing with long cable

IX-HLSH100
100W halogen lamp housing centerable

U-RMT
Extension cord

U-LH100-3
100W halogen lamp housing

IX-CI
Critical illuminator

U-FWR
Motorized reflected filter wheel

IX2-RFSS
Rectangular field stop

IX2-RFSPOT
Pinhole field stop

IX2-RFAL
L shape fluorescence illuminator

IX2-RFA
Fluorescence illuminator

IX2-RFAEVA*2
Total internal reflection fluorescence illuminator

32ND6-2
32ND12-2
32ND25-2
32ND50-2
32FR80
ND filters

U-DULHA
Dual lamp housing attachment

U-CLA
Collector lens adjustment

U-LH100HG
100W mercury lamp housing

U-LH100HGAPO
100W mercury apo lamp housing

Power supply unit

U-LH75XEAPO
75W xenon apo lamp housing

Power supply unit

FV5-FUR*2
Fiber unit

FV5-LAAR*2
Argon ion laser

IX2-PRBP1*2
100% prism unit for lower back port

IX2-PRBP8*2
20/80% prism unit for lower back port

IX2-PRS1*2
20/80% prism unit for side port

IX2-PRS8*2
100% prism unit for side port

IX2-LBPC*2
Lower back port unit with C mount unit

*1 Not available in some areas
*2 To consult with local Olympus subsidiary